

Histochemical Detection of GUS (beta-glucuronidase) (1)

This protocol is for the Histochemical Detection of GUS (beta-glucuronidase) (1)

Staining buffer, prepare immediately before use as follows:

Stock solutions	Volume per 1 ml staining buffer	Final concentration
Water, nuclease-free	830 µl	
1 M sodium phosphate (pH 7.0)	100 µl	0.1 M
0.5 M EDTA, pH 8.0	20 µl	10 mM
10% Triton X-100	10 µl	0.1% (v/v)
50 mM K ₃ Fe(CN) ₆	20 µl	1 mM
0.1 M X-Gluc (50 mg/ml) in dimethylformamide	20 µl	2 mM

Staining procedure for cells and tissues:

1. Remove media from the cells or tissue.
2. Immerse cells/tissue in fresh staining buffer.
3. Incubate cells/tissue for 12-24 hours at 37°C.
4. Remove the staining buffer.
5. Wash with several changes of 50% ethanol (up to 12 hours per wash), until the cells/tissue clears.
6. Count dark blue cells.

Reference

1. Jefferson, R., Assaying chimeric genes in plants: the GUS gene fusion system, *plant Mol Biol Rep*, 5, 387-405, 1987.

thermoscientific.com

© 2012 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

North America
Technical Services:
 techservice.genomics@thermofisher.com
Customer Services:
 customerservice.genomics@thermofisher.com
 Tel 800 235 9880
 Fax 800 292 6088

Europe and Asia
Technical Services:
 techservice.emea.genomics@thermofisher.com
Customer Services:
 customerservice.emea.genomics@thermofisher.com

Thermo
 S C I E N T I F I C

Part of Thermo Fisher Scientific