Introduction

The Varipette/Maxipettor® becomes a functional unit when a Varitip is attached. The Varitip consists of a piston, barrel and tip and holds the liquid to be dispensed.

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Design principle

(Fig.1)

- 1 Varipette/Maxipettor
 - 2 Varitip S/Maxitip S, dispensing part (barrel with piston) (2.50–10.0 ml)
 - 3 Maxitip
The Maxitip is attached to the dispensing part and is interchangeable.
- The Maxitip is autoclavable.
 - The Varitip S/Maxitip S with Maxitip is for pipetting in the range from 2.50 ml to 10.00 ml. The long Maxitip is designed for aspirating liquids from tall and long-necked vessels (diameter of tip = 10 mm, fits into flasks with standard ground joint 12/21). During pipetting, solution only enters the Maxitip, therefore, if the Maxitip is exchanged, samples and reagents can be dispensed in succession without carryover.

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Varitip / Maxitip Kit

Abb./Fig. 2 **Valve in Maxitip** (Fig. 2)

Can be inserted either way round into the Maxitip. The valve closes the Maxitip and prevents dripping.

The valve closes the Maxitip with its own weight (density = 1.78).

- Before filling, hold the Maxitip against the inside of the vessel so that the valve is pushed back once.
- During filling, the flowing liquid lifts the valve. When dispensing or moving the piston downwards, hold the Maxitip at an angle against the inside of the vessel so that the valve is pushed up.
- Hold the Varipette/Maxipettor with filled Maxitip and valve vertically.

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