

Electro Zap™ Solution

Store at room temperature.

Catalog #: AM9785

Product Description: A special formulation designed to eliminate RNase contamination from electrodes.

Amount: 250 mL Concentration: 1X strength

Storage Conditions: Store at room temperature.

USER INFORMATION

General Information: ElectroZap™ Electrode Decontamination Solution was formulated using three ingredients known to be active against

> RNases and ideal for treating electrodes. It effectively removes high levels of RNase contamination on contact with surfaces and leaves no residues that are inhibitory to enzymatic reactions. [Note: A different product, RNaseZap® Solution (Cat #AM9780), is available for treating non-electrode surfaces, including plastic and glassware. See

www.ambion.com for more information.]

The solution comes ready to use. Apply ElectroZap Solution by spraying it directly onto pH electrodes, thoroughly Applications:

coating the electrode's surface. Small parts may be cleaned by briefly soaking them in the solution. After

application, all treated surfaces should be rinsed thoroughly two times with distilled water.

Note: ElectroZap Solution is not compatible with corrodable metals (e.g. aluminum) or rubber (gaskets, etc.). It is

also not meant for treating solutions.

OTHER INFORMATION

Material Safety Data Sheets (MSDSs) can be printed or downloaded from product-specific links on our website at the following **Material Safety Data Sheets:**

address: www.ambion.com/techlib/msds. Alternatively, e-mail your request to MSDS_Inquiry_CCRM@appliedbiosystems.com. Specify the catalog or part number(s) of the product(s), and we will e-mail the associated MSDSs unless you specify a preference for fax delivery. For customers without access to the internet or fax, our technical service department can fulfill MSDS requests placed by

telephone or postal mail. (Requests for postal delivery require 1-2 weeks for processing.)

Warranty and Liability: For research use only. Not for use in diagnostic procedures.

Applied Biosystems is committed to delivering superior product quality and performance, supported by industry-leading global service

and technical support teams. Warranty information for the accompanying consumable product is available at

www.ambion.com/info/warranty in "Limited Warranty for Consumables," which is subject to the exclusions, conditions, exceptions, and limitations set forth under the caption "EXCLUSIONS, CONDITIONS, EXCEPTIONS, AND LIMITATIONS" in the full warranty statement. Please contact Applied Biosystems if you have any questions about our warranties or would like information about

Information in this document is subject to change without notice. Applied Biosystems assumes no responsibility for any errors that

may appear in this document

Applied Biosystems disclaims all warranties with respect to this document, expressed or implied, including but not limited to those of merchantability or fitness for a particular purpose. In no event shall Applied Biosystems be liable, whether in contract, tort, warranty, or under any statute or on any other basis for special, incidental, indirect, punitive, multiple or consequential damages in connection

with or arising from this document, including but not limited to the use thereof.

Trademarks, Patents, and

Licensing:

Applied Biosystems, AB (Design), and Ambion are registered trademarks, and Electro Zap is a trademark of Applera Corporation or its subsidiaries in the US and/or certain other countries. All other trademarks are the sole property of their respective owners.

© 2008 Ambion, Inc. All rights reserved. 4386651A